

Effects of washing and re-using autosampler vials on sample loss – Tech Information

As reported by ThermoFisher Scientific, laboratories may try to save money by washing their used autosampler vials and reusing them. However, these products are designed to be disposable and reuse can cause numerous QC issues, such as contaminant peaks, sample loss, and physical deformations on the glass surface. Here a variety of vials are tested for sample loss. Only the new vial showed zero loss for all samples. The others showed variable amounts of loss.



The variable nature of the sample loss can make identifying the cause of the issue problematic. You can avoid all this by simply using new vials each time.

Printed from the Chrom Resource Center

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 05-05-2024