



## Domoic Acid in Seafood Samples Analyzed by LCMS – AppNote

### Domoic Acid (DA) From Seafood Without Derivatization

Several Seafood and Plankton samples were tested for Domoic Acid (DA) which, as a highly polar compound, is poorly Retained on most Reversed Phase Columns. This developed Method based on HPLC and MS detection does not require derivatization of Domoic Acid.

The concentration of this potent toxin varied quite substantially between the samples (*from 40ng / mL in Plankton to 3700ng / mL in Mussels*). Despite different sample matrices, it was possible to detect and analyze Domoic Acid in a wide range of concentrations.

The results were very Repeatable (*see 5 overlaid injections in Figure F*). The detection limit depends on the instrument used. The Method can be used by regulatory agencies responsible for monitoring the occurrence of toxins such as the analysis of Domoic Acid in seafood samples.

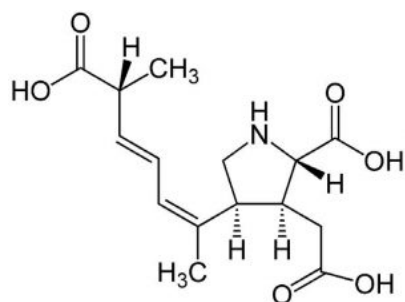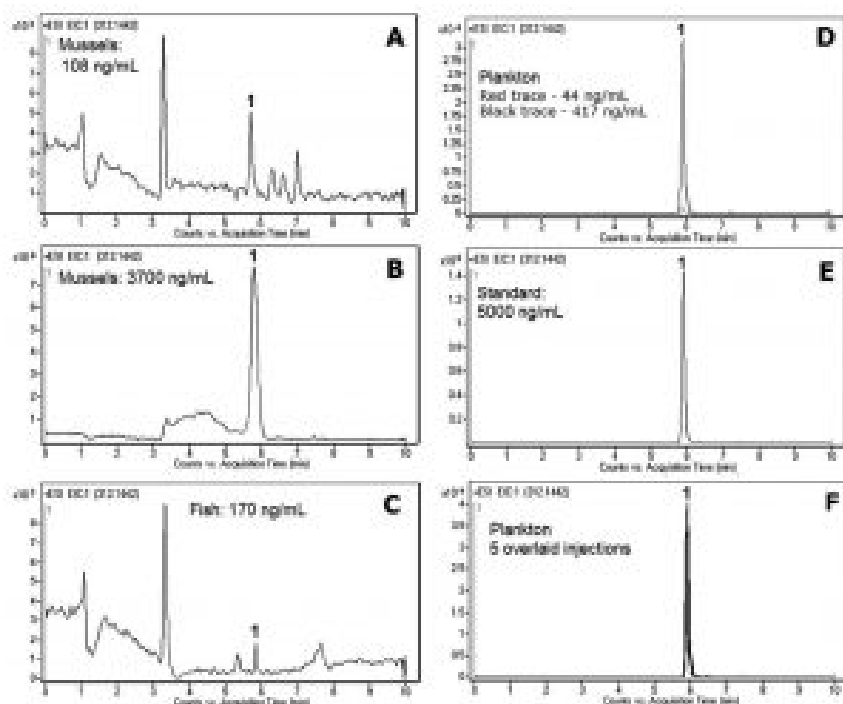

**Peak:**

Domoic Acid 312.1442 m/z (M+H)<sup>+</sup>

## Method Conditions

**Column:** Cogent Diamond Hydride™, 4µm, 100Å

**Catalog No.:** 70000-15P-2

**Dimensions:** 2.1 x 150mm

**Mobile Phase:**

A: 50% MeOH / 50% DI Water / 0.1% Formic Acid

B: Acetonitrile / 0.1% Formic Acid

**Gradient:**

Time (minutes)	%B
0	95
7	30
8	30
10	95

**Post Time:** 5 minutes

**Injection vol.:** 1µL

**Flow rate:** 0.4mL / minute

**Detection:** ESI - neg - Agilent 6210 MSD TOF Mass Spectrometer

**Sample Preparation:** Samples from Dr. R.Kudela, Univ. of California, Santa Cruz, Dept. of Marine Sciences.

Samples were injected as received.



## Attachment

**No 140 Domoic Acid in Seafood Samples Analyzed by LCMS pdf** 0.3 Mb [Download File](#)

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 05-21-2024