

## Is the Holmium Oxide Solution intended as a stand alone product or solely as a replacement solution for the HSQ Kit – FAQ

The Holmium Oxide solution (catalog number [PQA-HOX-01](#) or [PQA-HOX-A](#)) can be used as a stand-alone item without having to order the entire HSQ Kit. It is a component of the Kit but can be used to perform wavelength **accuracy** qualification on its own. You can also order it as a replacement solution if you already have the HSQ Kit™.

The solutions do not come with the Excel template included in the kit but it is not necessary for wavelength qualification. You simply pick any or all of the 14 bands from 241 – 641 nm and compare their found values to the official values provided on the supplied certificate of analysis.

For more detail, please refer to the below downloadable pdf instructions.



Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 05-18-2024