



## Why is it not recommended to use FEP tubing above 80°C when the melting point is much higher - FAQ

**FEP polymer has a melting point of 260°C and a general recommended maximum operating temperature during usage of around 200°C.**

If pressure is applied in the tubing, however, the maximum temperature should be lowered accordingly. The maximum temperature during usage should be 80°C under these conditions. If no pressure is applied, the tubing can be used up to 200°C.

Click [HERE](#) for low pressure HPLC Tubing ordering information and pictures.

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 05-05-2024