


## Cogent Diamond Hydride 2.0 HPLC column retention time precision – AppNote

This graph is from a study investigating the day to day retention time precision with two test solutes on the Cogent Diamond Hydride 2.0™ HPLC column

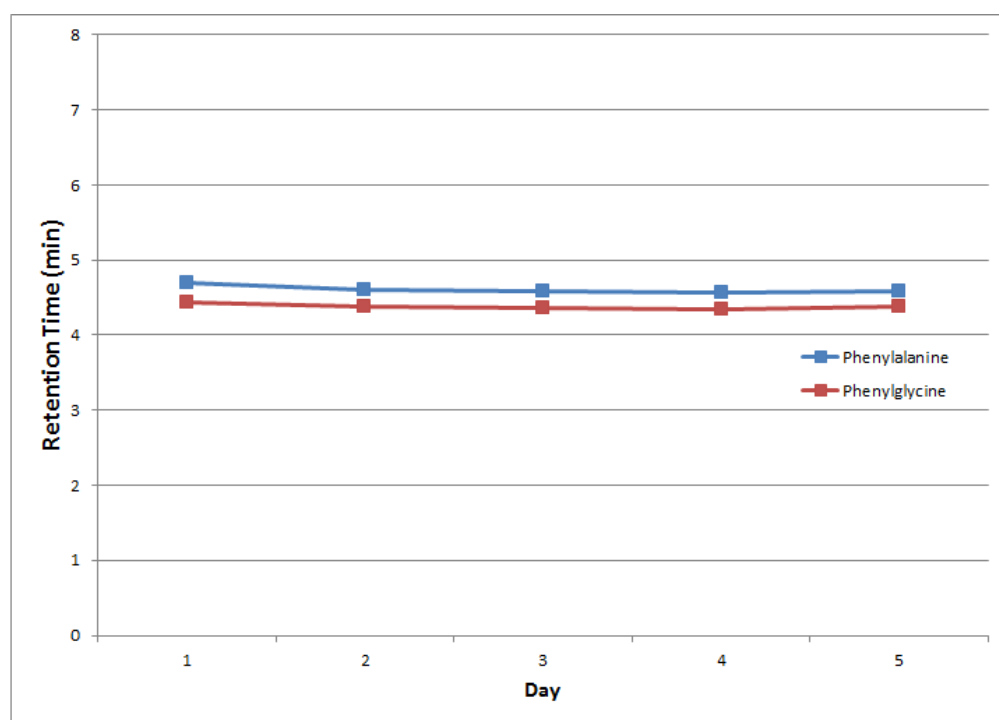

### Peaks:

Blue Trace: Phenylalanine

Red Trace: Phenylglycine

Many factors can affect column **robustness** and lifetime, so it is not possible to define specific values for those parameters.

### Mobile Phase:

A: DI Water / 0.1% Formic Acid

B: Acetonitrile / 0.1% Formic Acid

**Flow Rate:** 0.3mL / minute

**Injection Volume:** 0.2uL

**Detection:** UV 254nm

### Gradient:

Time (minutes)	%B
0	90
6	20
7	20



## [Cogent Diamond Hydride HPLC Column Ordering Information](#)

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 05-20-2024