


Atorvastatin drug product analyzed with HPLC – AppNote


API well separated from degradation product

Using this Simple Linear Gradient, Atorvastatin is well separated from its main degradation Product. The degradant is formed under acidic conditions and therefore its Separation is important for Stability Indicating Methods for pharmaceutical formulations of Atorvastatin. For Methods requiring non-degrading conditions, it is important to choose diluent and Mobile Phase conditions with pH sensitivity in mind.

As such, Ammonium Acetate was chosen for the diluent and Mobile Phase additives in order to avoid any degradation from these sources. Data obtained under non-degrading conditions is shown in *Figure A*. An overlay of five consecutive runs of the degraded extract is shown in *Figure B*. The Repeatability of the Method is demonstrated by the consistent retention time (%RSD of 0.10).

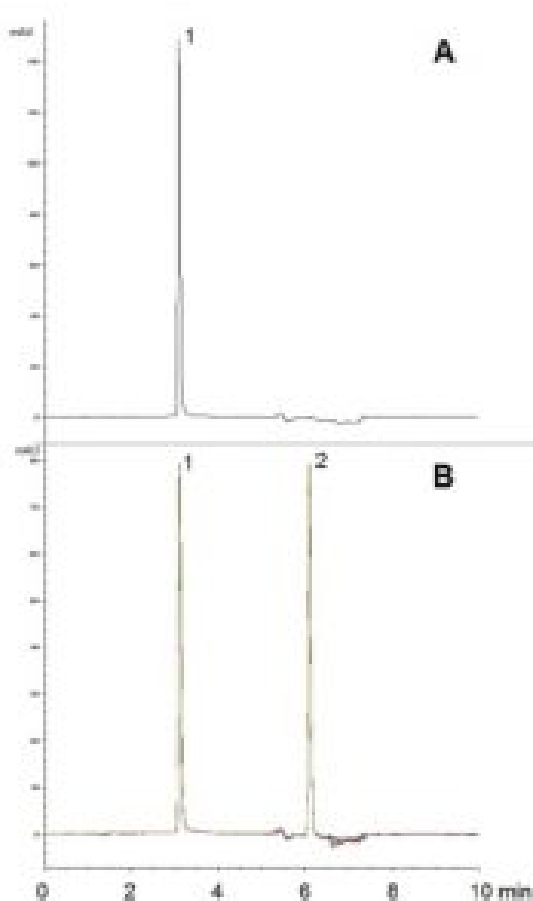

1. Atorvastatin
2. Degradant

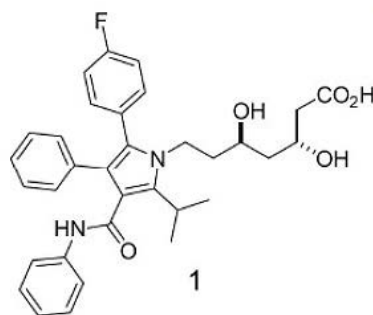

Peaks:

Method Conditions

Column: Cogent Bidentate C18™, 4µm, 100Å

Catalog No.: 40018-15P

Dimensions: 4.6 x 150mm

Printed from the Chrom Resource Center
Copyright 2024, All Rights Apply

MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451
tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Mobile Phase:

A: DI Water / 10mM Ammonium Acetate

B: 90% Acetonitrile / 10% DI Water / 10mM Ammonium Acetate

Gradient:

Time (minutes)	%B
0	40
5	100
6	40

Injection vol.: 2µL

Flow rate: 0.8mL / minute

Detection: UV @ 248nm

Sample Preparation:

A: Atorvastatin tablet extract

B: Degraded Atorvastatin tablet extract

40mg strength tablet ground was placed in 100mL volumetric flask. 50mL of Mobile phase B was added to flask. Sample was vortexed for 5 minutes, sonicated for 5 minutes, then diluted to mark with Solution A. A portion was filtered through a 0.45µm Nylon Syringe Filter (MicroSolv Tech Corp.). Final Working Solution was then diluted 4x in **Sample A)** 50:50 A / B or **Sample B)** 1.0 N HCL in 50:50 A / B.

Note: Atorvastatin is a competitive inhibitor of 3-Hydroxy-3-Methylglutaryl Co-enzyme A (HMGCoA) reductase, which catalyzes the rate-limiting step in Cholesterol biosynthesis. As such, Atorvastatin is used to reduce plasma levels of low-density lipoprotein (LDL) Cholesterol, which are known to contribute to the development of Atherosclerosis.

**Attachment**

Atorvastatin Drug Product Analyzed with HPLC pdf [Download File](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com