
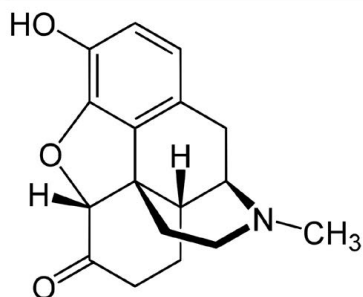


## Hydromorphone Tablet Analyzed with HPLC – AppNote

### Alternative to USP Method for this Opioid

The USP Method for Hydromorphone Tablet Assay calls for the use of Sodium Dodecyl Sulfate (SDS) in the Mobile Phase. In addition to not being LCMS compatible, adding it is time consuming to load and unload from the HPLC column. Often the columns are permanently altered and cannot be used for other analyses.

The Method conditions presented here do not require SDS and can be used in LCMS, which is advantageous since Hydromorphone is often analyzed in blood or urine matrices. Data from two column lots is shown to illustrate Method Robustness and Consistency.

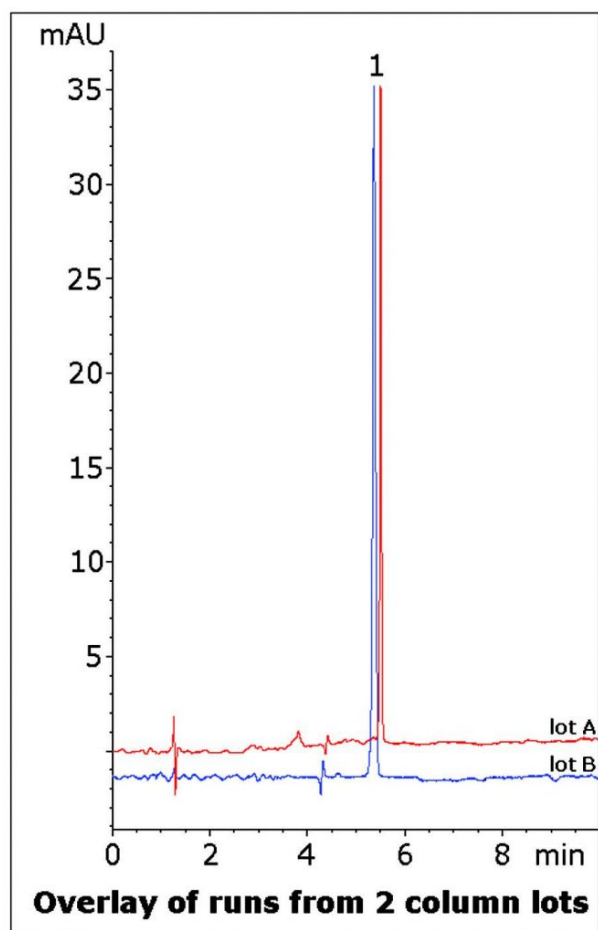

Peak:

Printed from the Chrom Resource Center  
Copyright 2024, All Rights Apply  
**MicroSolv Technology Corporation**  
9158 Industrial Blvd. NE, Leland, NC 28451  
tel. (732) 380-8900, fax (910) 769-9435  
Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)  
Website: [www.mtc-usa.com](http://www.mtc-usa.com)

## Method Conditions

**Column:** Cogent Diamond Hydride™, 4µm, 100Å

**Catalog No.:** 70000-7.5P

**Dimensions:** 4.6 x 75mm

**Mobile Phase:**

A: DI Water with 0.1% Formic Acid (v/v)

B: Acetonitrile with 0.1% Formic Acid (v/v)

**Gradient:**

Time (minutes)	%B
0	95
1	95
6	40
7	95

**Post Time:** 3 minutes

**Injection vol.:** 2µL

**Flow rate:** 1.0mL / minute

**Detection:** UV @ 280nm

**Sample Preparation:** 2mg strength Hydromorphone Tablet was ground and added to a 10mL volumetric flask. A portion of 50:50 Solvent A / Solvent B diluent was added and the flask was sonicated 10 minutes. It was then diluted to mark and mixed. A portion was filtered with a 0.45µm Nylon Syringe Filter (MicroSolv Tech Corp.).

**t<sub>0</sub>:** 0.9 minutes

**Note:** Hydromorphone is a derivative of Morphine and is used as an analgesic to treat moderate to severe pain. It is available by prescription under various trade names such as Palladone® and Dilaudid®.



### Attachment

No 248 Hydromorphone Tablet Analyzed with HPLC pdf 0.4 Mb [Download File](#)