
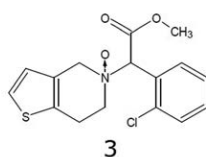
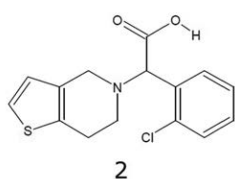
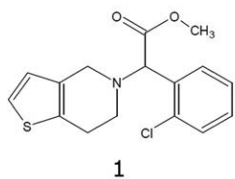
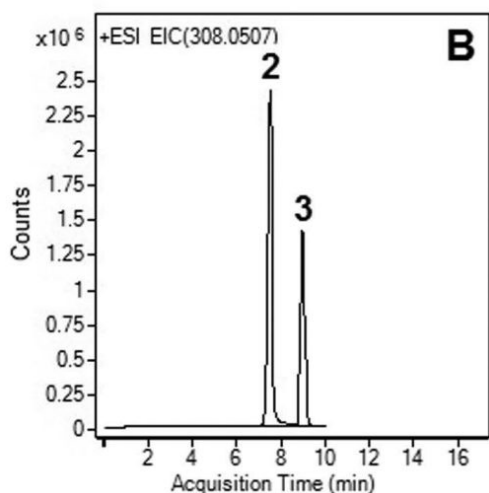


## Analysis of Clopidogrel with LCMS – AppNote

### Analysis of the API and Degradants

A Sensitive, Selective, and rapid LCMS Method was developed for the simultaneous Quantification of Clopidogrel (Plavix®) and degradants.

This Method can be used for the determination of Clopidogrel in commercial tablets for Quality Control with an application for content uniformity test. The Method is also stability-indicating as it is suitable for the determination of the API in the presence of its degradation products under all stress conditions using HCL, NaOH, light and Hydrogen Peroxide.

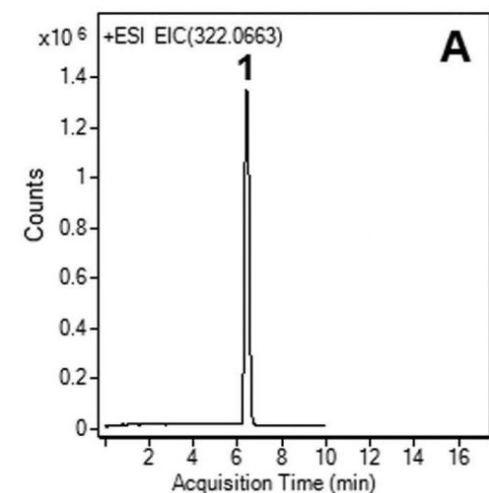

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

#### Peaks:

1. API: Clopidogrel, m/z 322.0663 [M+H]<sup>+</sup>

2. Degradant: Clopidogrel Acid, m/z 308.0507 [M+H]<sup>+</sup>

3. Degradant: Clopidogrel N-oxide, m/z 338.02 [M+H]<sup>+</sup>

## Method Conditions

**Column:** Cogent Diamond Hydride™, 4µm, 100Å

**Catalog No.:** 70000-15P-2

**Dimensions:** 2.1 x 150mm

### Mobile Phase:

A: DI Water / 0.1% Formic Acid (v/v)

B: Acetonitrile / 0.1% Formic Acid (v/v)

### Gradient:

Time (minutes)	%B
0	95
2	95
7	60
8	95

**Temperature:** 25°C

**Post Time:** 3 minutes

**Injection vol.:** 1µL

**Flow rate:** 0.4mL / minute

**Detection:** ESI - POS - Agilent 6210 MSD TOF Mass Spectrometer

**Sample Preparation:** 50mg strength Plavix® tablet was ground and diluted in 50:50 Solvent A / Solvent B mixture to 50mL. The solution was sonicated and filtered through a 0.45µm Nylon Syringe Filter (MicroSolv Tech Corp.).

**t<sub>0</sub>:** 0.9 minutes

### Figures:

*Fig. A: Non-degraded extract: The stock solution was diluted 1:10 with 50:50 Solvent A / Solvent B mixture. Only the API peak is observed.*

*Fig. B: Base degradation with heating: The stock solution was diluted 1:10 with 50:50 1N NaOH / Acetonitrile mixture and then heated at 85°C for 30 minutes. The API peak is no longer observed but degradants (Peak 2 and 3) are now present.*



### Attachment

No 258 Forced Degradation of Clopidogrel Analysis LC-MS pdf 0.3 Mb

[Download File](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)