



Using Buffer Additives in CZE

When you add Modifiers (Buffer Additives) to your CZE method, the following changes can be expected:

Joule Heating: Adding high concentrations of organic modifiers or other additives can cause an increase in current or an increase in Joule Heating.

Viscosity: Adding high concentrations of additives such as methylcellulose will increase viscosity. Also this can cause a molecular sieving effect as well.

Electro Osmotic Flow: Depending on the additive, the EOF can be suppressed or enhanced but organic solvents usually lower the EOF.

Current: When using high concentrations of surfactants as an additive, this can cause a high current to be produced.

Analyte to Wall Interaction: Depending on the additive, this can prevent analyte to wall interaction or binding to the wall.

Migration Time: Depending on the additive, you can increase or decrease migration times. Additives are also used to change migration order.

Resolution: Additives can enhance your resolution and change your peak shape. Using additives permit indirection detection to take place.

Electrophoretic Mobility: Using additives can enhance analyte electrophoretic mobility.

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com