
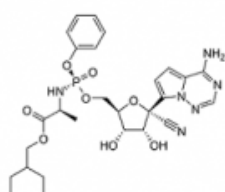


## Remdesivir and GS-441524 a ProDrug – AppNote


### Retention of polar and non-polar drug compounds in ANP

Remdesivir, an antiviral drug, and GS-441524, a pro-drug converted in the body as a ribonucleotide analog, is as a mixture that may be difficult to analyze in routine chromatographic analysis. As GS-441524 is polar, it makes it more difficult to retain in typical Reversed Phase conditions.

By use of the Cogent Diamond Hydride™ Column in ANP, we have illustrated how this contrasting mixture of compounds can be retained and still possess excellent peak shape. This method demonstrates an easy HPLC method that can be easily transferred to LCMS.

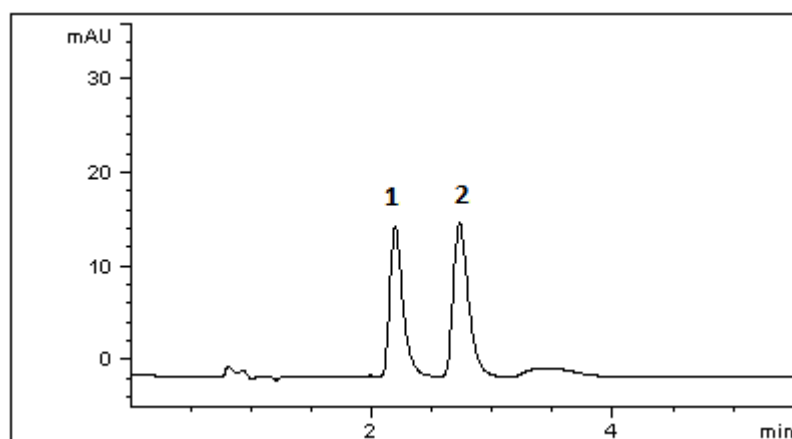

Remdesivir

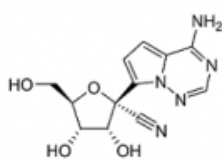

GS-441524

### Peaks:

1. Remdesivir
2. GS-441524

### Method Conditions:

**Column:** Cogent Diamond Hydride™, 4 μm, 100 Å

**Catalog No.:** 70000-75P

**Dimensions:** 4.6 x 75mm

### Mobile Phase:

- A. DI Water / 0.1% Formic Acid (v/v)
- B. Acetonitrile / 0.1% Formic Acid (v/v)

**Isocratic:** 92% B

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

**Injection Volume:** 2ul

**Flow Rate:** 1.0 mL / minute

**Detection:** 254 nm

**Samples:** Remdesivir and GS-441524 in 0.1 mg/mL concentration

**Note:** Remdesivir is a broad-spectrum antiviral drug that is able to diffuse into cells where it is converted to GS-441524 mono-phosphate. This is achieved via the actions of esterases and a phosphoamidase; which in turn is further phosphorylated to its active metabolite triphosphate by nucleoside-phosphate kinases GS-441524. GS-441524 has previously been researched as a treatment for feline infectious peritonitis (FIP), a lethal coronavirus disease which affects domestic cats.



## Attachment

**Remdesivir and GS-441524 App Note.pdf** 0.1 Mb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)