



## Sample diluents should match your starting mobile phase – Tips & Suggestions

For best results, it is always best to try to dissolve your samples in as close as possible to your isocratic **eluent** or the start of your chosen gradient.

If this is not possible gradually increase the percentage of the solvent in your **eluent** in which the target is most soluble.



Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)