



## Injection volume suggestions for Cogent TYPE-C HPLC columns - Tips & Suggestions

When using an analytical Cogent TYPE-C™ column it is recommended to inject 1uL to 10uL of the sample which is prepared either in a diluent that matches, as closely as possible, the starting mobile phase composition of a gradient method or the method's isocratic mobile phase.

Larger volume injections that are performed on the columns are probably overloading the columns.



Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 07-03-2024