



## Storing the HPLC column supplied in the HSQ kit – How To

One of the best features of the MicroSolv HSQ Kit™ is that the kit was designed for ease and speed. The certified, special HPLC column is shipped and stored in the same mobile phase that is used during the testing.

The column is always ready to go with minimal equilibration time. The mobile phase / storage solvents were chosen for long column life. The solutions should be stored at **room temperature** in a dry place and never refrigerated.

Click [HERE](#) for PQ Kit™ and HSQ Kit™ ordering information and pictures.

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)