



## What is the mobile phase used in the HSQ Kit and how much is used - FAQ

**For isocratic system testing, prepare one liter of 13:87 acetonitrile:0.1% acetic acid (1mL/L).**

For gradient testing, prepare 2 liters.

Click [HERE](#) for PQ Kit™ and HSQ Kit™ ordering information and pictures.

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 07-22-2024