



## Can I autoclave the MicroSolv AQ brand 25mm PTFE syringe filters - FAQ

**Yes, under conditions of 121°C for a maximum of 15minutes, the PTFE AQ™ brand of syringe filters can be autoclaved.** Maximum operating temperature (during continuous continuous autoclaving) we suggest that you do not exceed 80°C to maintain optimum performance.

Click [HERE](#) for ordering information and pictures of AQ™ brand of PTFE syringe filters

Attachment:

MicroSolv Filters Equivalency Study pdf [Download File](#)

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 07-22-2024