

What determines the retention of a molecule in reversed phase HPLC – FAQ

There has been much research and work done in this area and many books have been written about this topic.

However, to give a very short answer you could say that much of the retention of a column in **Reversed Phase RP** HPLC is due to what is commonly called hydrophobicity. This natural phenomena is due mostly to the size of the hydrophobic (water resistant) area of a molecule. Retention will increase with the amount of water in the mobile phase. In general, the more hydrophobic the molecule the longer it should be retained.

If a molecule is “solvo-phobic” the molecule will retain elsewhere, namely the column. This goes to the old adage in chromatography “like likes like”.

NOTE: This does not apply to Normal Phase HPLC, Aqueous Normal Phase ANP or Ion Exchange Chromatography.



Printed from the Chrom Resource Center

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 04-05-2024