



## Analyzing metformin in plasma samples how do I obtain reproducible results and peak shapes - FAQ

**Suggestion:** Try substituting formic acid in DI water and adding ammonium acetate (*10mM*) and acetic acid (*1%*) to the aqueous phase. This should improve the reproducibility of peak areas and retention times.

**In addition,** extensive reconditioning of the analytical column with pure acetonitrile in between runs is essential for good reproducibility when working with plasma samples.



Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 04-05-2024