
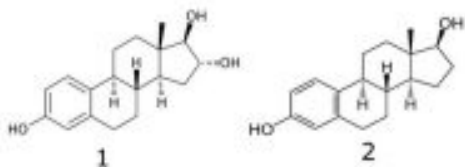


Separation of Estriol, Estradiol, & Progesterone

This gradient method features a separation of the three components of a hormone replacement formulation. *Figure A* shows a five run overlay of the formulation extract injections. *Figure B* shows a zoomed-in view so that the Estriol and Estradiol Peaks, which are present in much lower concentration than Progesterone, can be seen clearly. *Figure B* also shows separation of an impurity from the Progesterone Peak.

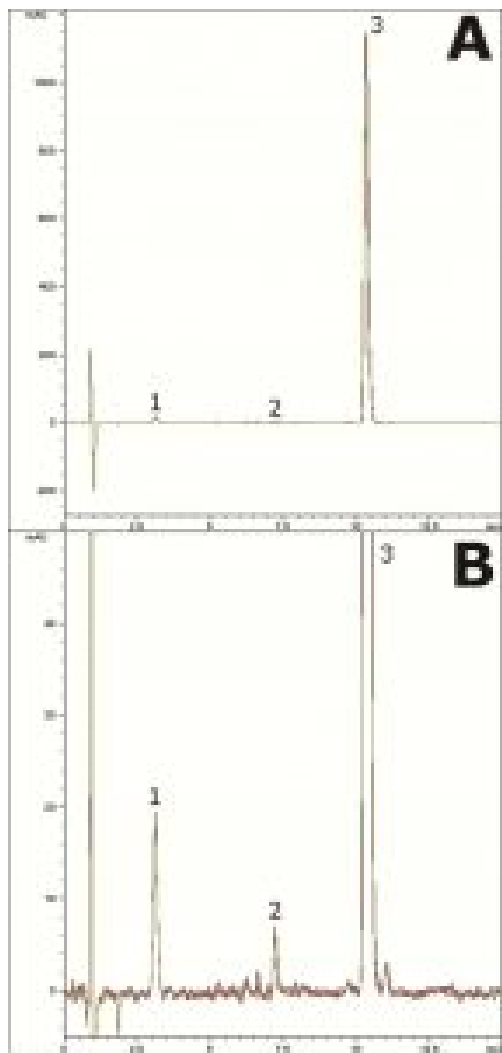

Peak:

1. Estriol
2. Estradiol
3. Progesterone

Method Conditions

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Column: Cogent UDC Cholesterol™, 4µm, 100Å

Catalog No.: 69069-7.5P

Dimensions: 4.6 x 75mm

Mobile Phase:

A: DI Water / 0.1% Formic Acid (v/v)

B: Acetonitrile / 0.1% Formic Acid (v/v)

Gradient:

Time (minutes)	%B
0	33
2	33
11	65
12	33

Post Time: 3 minutes

Flow rate: 1.0 mL / minute

Detection: UV @ 210nm

Injection vol.: 1µL

Sample Preparation: The contents of a capsule containing 0.124 mg Estradiol, 1.001 mg Estriol, and 50 mg Progesterone were added to a 25 mL volumetric flask. The flask was diluted to mark with solvent B and sonicated 10 minutes. Then A portion was filtered with a 0.45µm Nylon Syringe Filter (MicroSolv Tech. Corp.). Peak identities were confirmed by individual standards of 0.1 mg / mL in a Solvent B diluent.



Attachment

No 175 Hormone Replacement Capsule Analysis pdf 0.4 Mb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com