



What is the best way to regain flow in the capillary if I allowed it to dry out overnight - FAQ

The best way to regain flow is to remove 1 or 2mm of each end of the capillary. Then gently force methanol through the capillary using a syringe pump.

Click [here](#) to view a convenient hand pump for this purpose.

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 09-21-2024