


How do I know if a crimped cap on a vial has a proper seal – FAQ

The best way to know if you have a sufficient seal to prevent evaporation or leaking is to take the crimped vial and tightly hold it in your left hand. With your right hand, try to spin the crimped cap on the vial. If it moves even a little, you should remove the cap and adjust the crimping tool as shown below.

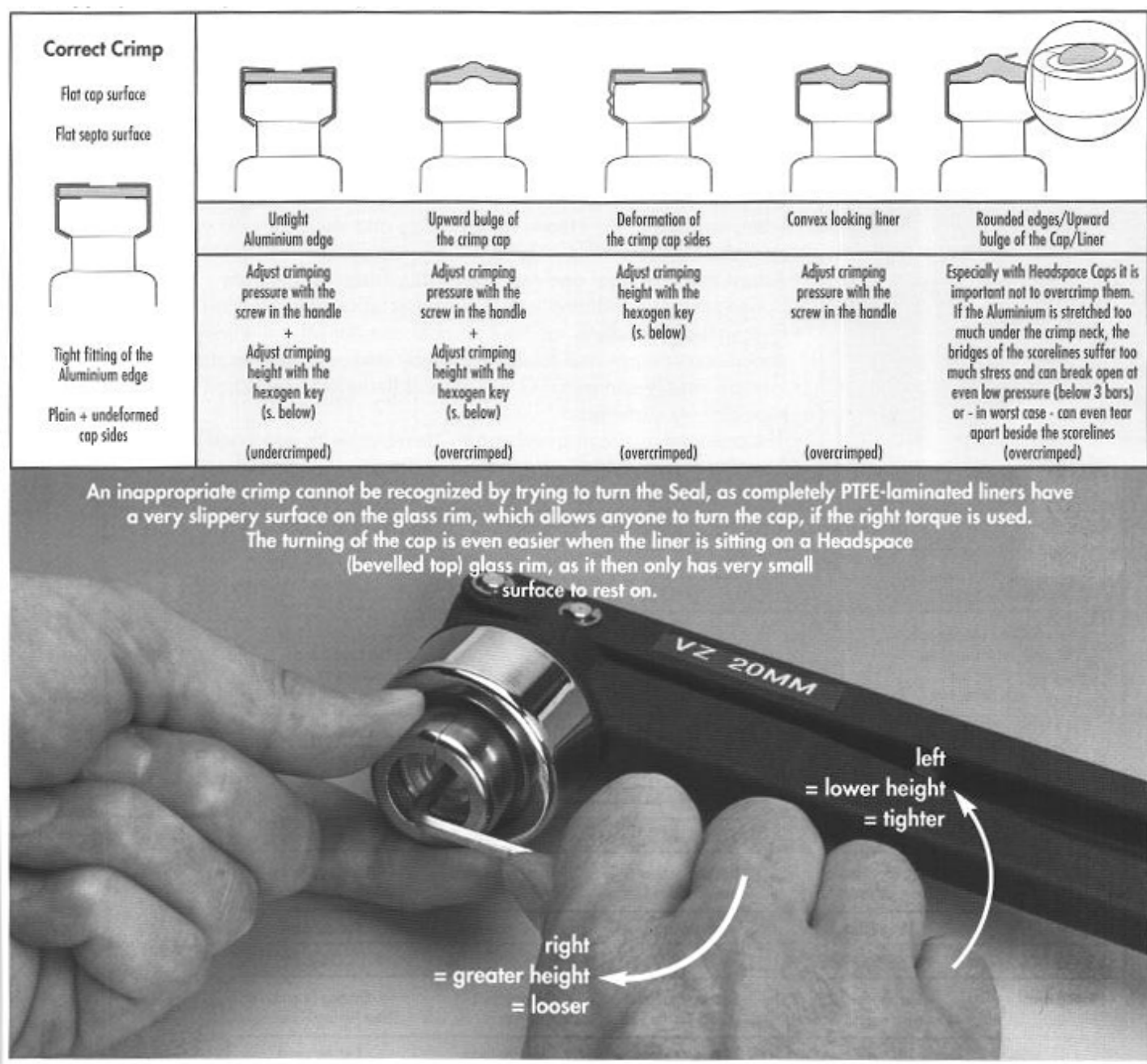

Printed from the Chrom Resource Center

Click [HERE](#) for Crimping Tool ordering information and pictures

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 05-19-2024