


EDTA Analysis with HPLC – AppNote

EDTA does not have a significant Chromophore, so to achieve UV Detection, in the Method shown below we used a pre-Column reaction of a Solution of Ferric Chloride with the Sample. The resulting EDTA/Fe³⁺ has significant UV Absorbance making this a very Sensitive Method.

Ethylenediaminetetraacetic acid is extremely difficult to analyze by itself however in its complexed form, it chromatographs well from matrices such as river sediment and other solutions.

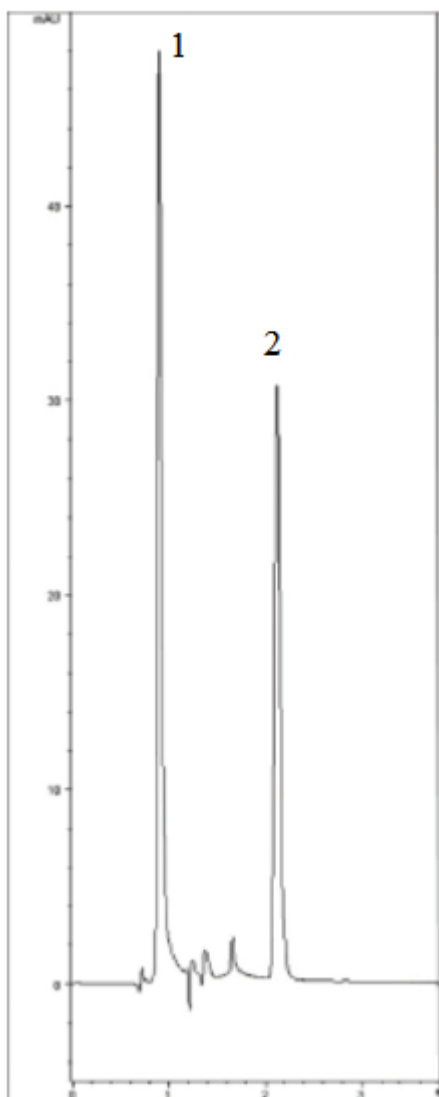


Peaks:

1. Water (solvent front)
2. EDTA Fe³⁺

Printed from the Chrom Resource Center
Copyright 2024, All Rights Apply

MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451
tel. (732) 380-8900, fax (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com

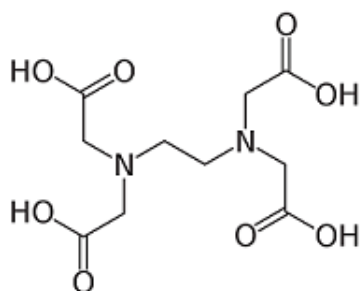

Ethylenediaminetetraacetic acid

Method Conditions:

Column: Cogent HPS C8™, 5µm, 120Å

Catalog No.: 75008-15P

Dimensions: 4.6 x 150mm

Mobile Phase: 98:2 DI H₂O/ Acetonitrile with 0.1% Acetic Acid (pH 3.5/2g/L Tetrabutylammonium Sulfate)

Temperature: 40°C

LOQ: 0.2µg / mL

Injection vol.: 20µL

Flow rate: 2mL / minute

Note: EDTA is a synthetic metal complexing reagent that is used in a wide variety of industrial applications. Used as a preservative, it has very low biodegradability thus remains in the environment for long periods of time. Found in sewer water, freshwater and ground water, it re-solubilizes precipitated toxic metals back into solution where they can be ingested by plants and animals.



Attachment

A74. EDTA Analysis with HPLC pdf 8.7 Kb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com