



## Digoxin Analyzed with HPLC - AppNote

### Forced Degradation Method

Digoxin had several Sugar Residues that can be hydrolyzed in acidic conditions. In this Forced Degradation Method, six degradants are observed. Some of these Peaks are likely the various hydrolyzed forms, including the Aglycone Digoxigenin.

The Method is capable of Separating a wide variety of Degradants. Figure A shows the Chromatogram of the non-degraded sample while Figure B shows 5 consecutive runs of the acid degraded extract.

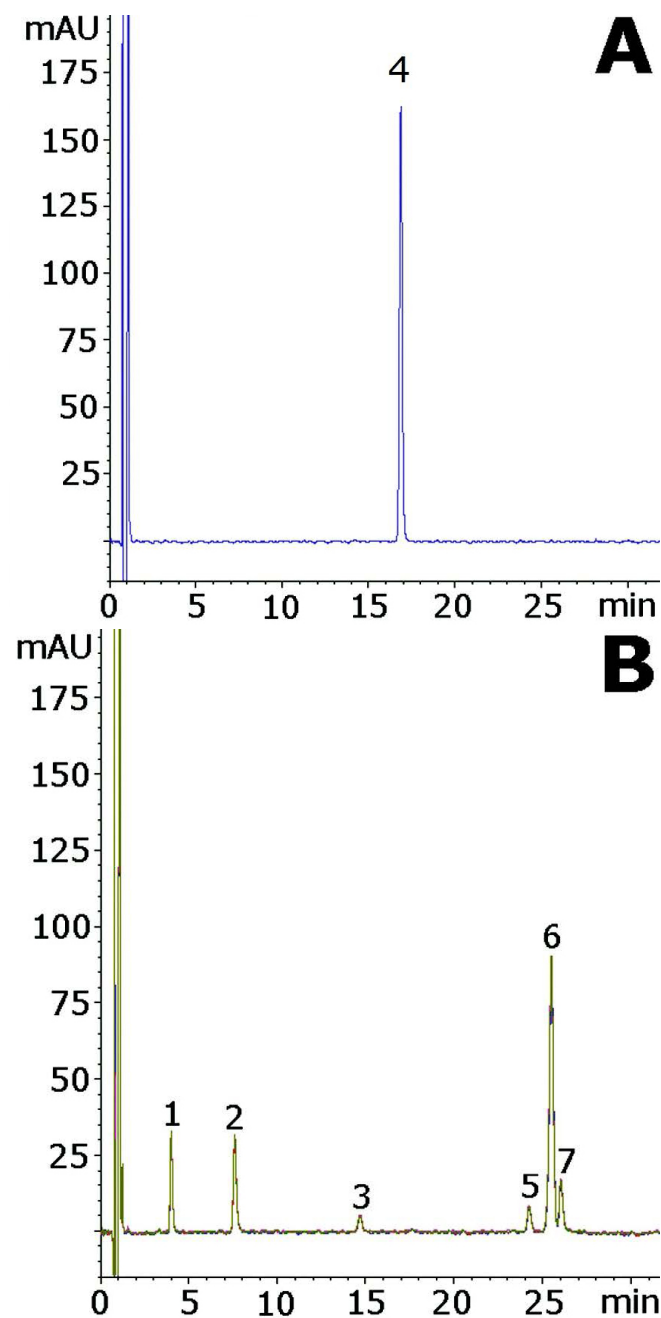

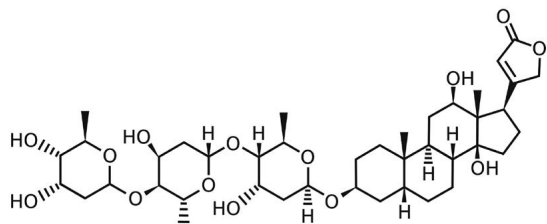

## Peaks:

1-3. Degradants

4. Digoxin

5-7. Degradants

## Method Conditions

**Column:** Cogent Bidentate C18™, 4µm, 100Å

**Catalog No.:** 40018-75P

**Dimensions:** 4.6 x 75mm

### Mobile Phase:

A: DI Water / 0.1% Formic Acid (v/v)

B: Acetonitrile / 0.1% Formic Acid (v/v)

### Gradient:

Time (minutes)	%B
0	15
2	15
26	37
28	80
29	15

**Post Time:** 3 minutes

**Injection vol.:** 10µL

**Flow rate:** 1.0mL / minute

**Detection:** UV @ 218nm

### Sample Preparation:

*Figure A:* Non-Degraded - 0.1mg / mL Digoxin in Methanol Diluent.

*Figure B:* Acid Degradation - 0.1mg / mL Digoxin in 50:50 Methanol / 1N HCL Diluent. Sample was heated at 85°C for 10 minutes.

**t<sub>0</sub>:** 0.9 minutes

**Note:** Digoxin is a cardiac glycoside used to treat a number of heart conditions. It is a natural product obtained from the Foxglove Plant (*Digitalis Purpurea*). Accurate Quantitation of Digoxin is crucial since it is very toxic at higher levels.



**Attachment**

**No 177 Digoxin Analyzed with HPLC pdf** 0.8 Mb [Download File](#)

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 07-22-2024