



Domoic Acid in Seafood Samples Analyzed by LCMS – AppNote

Domoic Acid (DA) From Seafood Without Derivatization

Several Seafood and Plankton samples were tested for Domoic Acid (DA) which, as a highly polar compound, is poorly Retained on most Reversed Phase Columns. This developed Method based on HPLC and MS detection does not require derivatization of Domoic Acid.

The concentration of this potent toxin varied quite substantially between the samples (*from 40ng / mL in Plankton to 3700ng / mL in Mussels*). Despite different sample matrices, it was possible to detect and analyze Domoic Acid in a wide range of concentrations.

The results were very Repeatable (*see 5 overlaid injections in Figure F*). The detection limit depends on the instrument used. The Method can be used by regulatory agencies responsible for monitoring the occurrence of toxins such as the analysis of Domoic Acid in seafood samples.

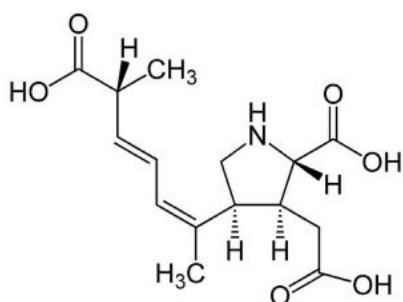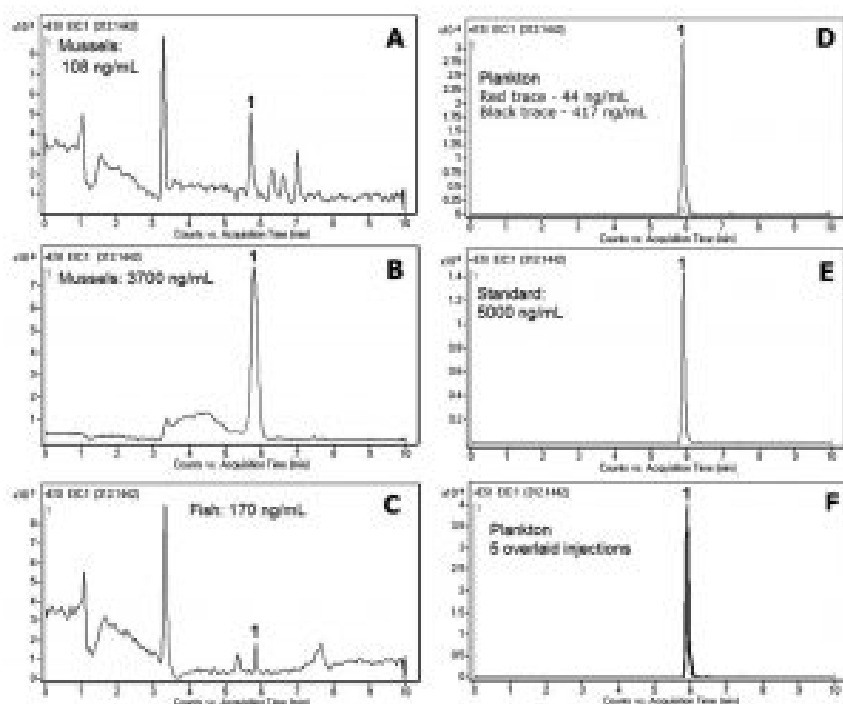

Peak:

Domoic Acid 312.1442 m/z (M+H)⁺

Method Conditions

Column: Cogent Diamond Hydride™, 4µm, 100Å

Catalog No.: 70000-15P-2

Dimensions: 2.1 x 150mm

Mobile Phase:

A: 50% MeOH / 50% DI Water / 0.1% Formic Acid

B: Acetonitrile / 0.1% Formic Acid

Gradient:

Time (minutes)	%B
0	95
7	30
8	30
10	95

Post Time: 5 minutes

Injection vol.: 1µL

Flow rate: 0.4mL / minute

Detection: ESI - neg - Agilent 6210 MSD TOF Mass Spectrometer

Sample Preparation: Samples from Dr. R.Kudela, Univ. of California, Santa Cruz, Dept. of Marine Sciences.
Samples were injected as received.



Attachment

No 140 Domoic Acid in Seafood Samples Analyzed by LCMS pdf 0.3 Mb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com