
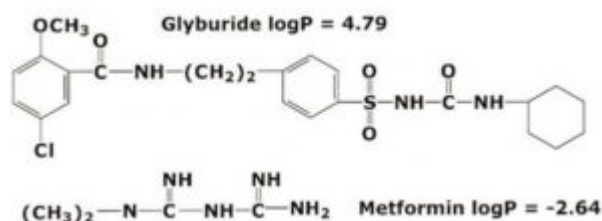
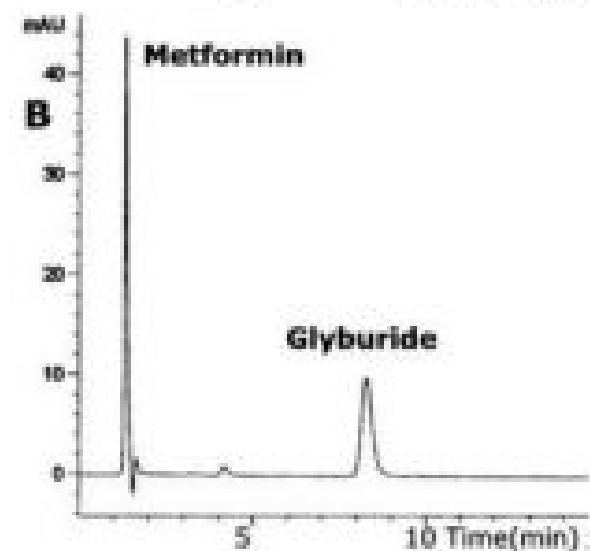


Glyburide and Metformin Analyzed with HPLC - AppNote

Separation of "Highly Polar" and "Non-Polar" Compounds in one Isocratic run

In this Method, the polar compound, Metformin, and the nonpolar compound Glyburide, can be retained on the same Stationary Phase (see A & B). Depending on the Mobile Phase composition either Metformin or Glyburide can be retained longer.

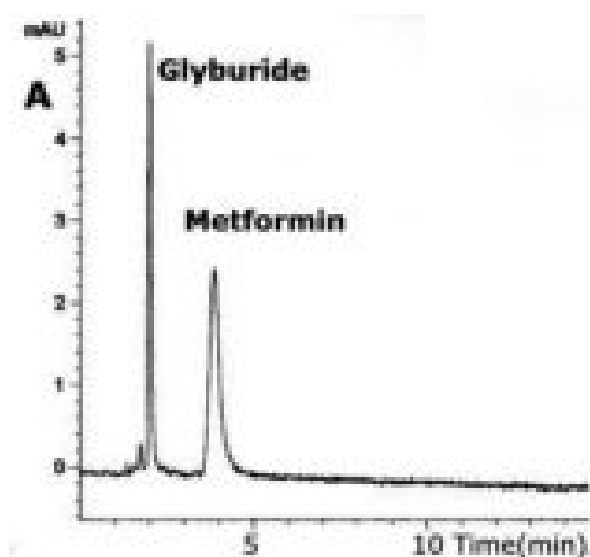

Method Conditions

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

MICROSOLV

Column: Cogent Bidentate C18™, 4µm, 100Å

Catalog No.: 40018-75P

Dimensions: 4.6 x 75mm

Mobile Phase:

A: 15% DI Water / 85% Acetonitrile / 0.5% Formic Acid

B: 50% DI Water / 50% Acetonitrile / 0.5% Formic Acid

Injection vol.: 1µL

Flow rate: 0.5mL / minute

Detection: UV @ 254nm

Sample Preparation: Stock Solution: 100µg / µL Glyburide and Metformin

Notes: Elution order was confirmed by LCMS, APCI+, with single ion monitoring Metformin (m/z 130) and Glyburide (m/z 369).



Attachment

No 08 Glyburide and Metformin Analyzed with HPLC pdf 0.2 Mb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com