
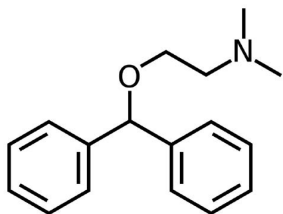


## Diphenhydramine HCL Capsule Analyzed with HPLC – AppNote

### Assay Method for Diphenhydramine HCL without Ion Pairing Agents

Diphenhydramine has a tertiary amine functional group that can produce tailing with conventional Columns in Reversed Phase HPLC. The USP Method features a Triethylamine additive in the Mobile Phase for this reason. These additives often take a significant time to completely load onto the Column and therefore adversely affect throughput and Robustness.

The Method in this Application Note produces an excellent Peak Shape using only Ammonium Acetate as a Mobile Phase Additive. In addition, the Reproducibility and Robustness of this Method is demonstrated through the overlay of four Chromatograms with Columns from four different batches.

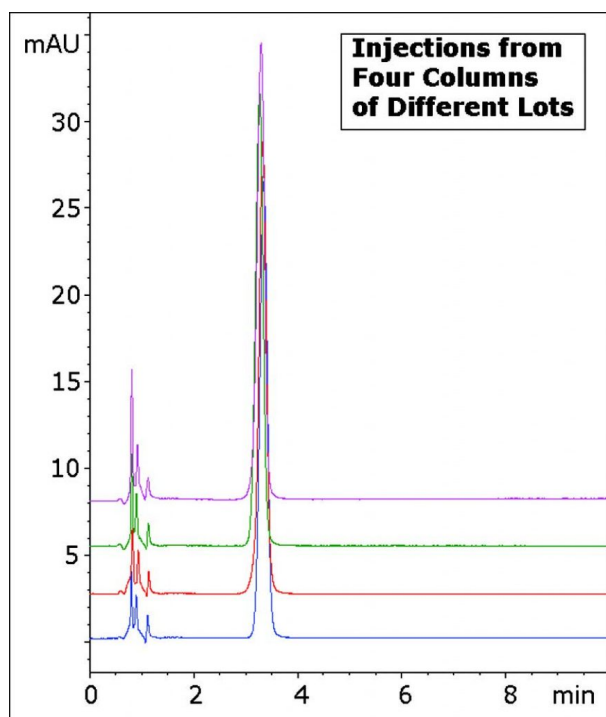

#### Peak:

Diphenhydramine

### Method Conditions

**Column:** Cogent Silica-C™, 4μm, 100Å

**Catalog No.:** 40000-7.5P

**Dimensions:** 4.6 x 75mm

**Mobile Phase:** 50:50 DI Water / Acetonitrile with 5mM Ammonium Acetate

**Injection vol.:** 2µL

**Flow rate:** 1.0mL / minute

**Detection:** UV @ 254nm

**Sample Preparation:** 25mg strength Benadryl® (Diphenhydramine) capsule was opened and placed in a 10mL volumetric flask with a portion of the Mobile Phase as diluent. It was sonicated 10 minutes and diluted to mark. Then a portion was filtered with a 0.45µm Nylon Syringe Filter (MicroSolv Tech Corp.).

**t<sub>0</sub>:** 0.9 minutes

**Note:** Diphenhydramine is a first generation antihistamine used primarily to treat allergies. It also has a significant sedative property, which is sometimes an undesirable side effect of its intended use. However, it is used in many formulations as a sleep aid as well.



## Attachment

**No 194 Diphenhydramine HCL Capsule Analyzed with HPLC.pdf** 0.4 Mb [Download File](#)

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 03-05-2024