


## What are the thread specifications for the PEEK columns - FAQ

The PEEK column hardware has *two specifications* for threading:

1. The female outlet threading found on the outside of the end fittings, where the end plugs or extra-column tubing ferrules connect. This threading is standard **10-32** for all items regardless of internal diameter (ID). This accommodates capillary tubing that is 1/16" outer diameter (OD).
2. The threading that connects the column tube to the end fittings. This varies depending on ID and is **1/4 - 28** for 2mm ID columns, **1/2 - 20** for 4.6mm, and **5/8 - 18** for 7.5mm.

### PEEK Hardware Ordering Information

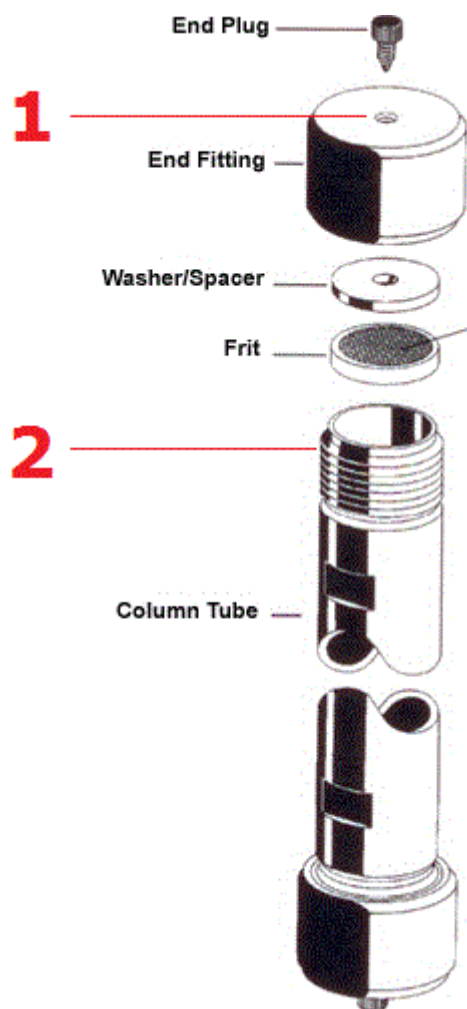

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 07-22-2024