

Sparging stones should not be touched directly by bare human hands if using with HPLC.

Always wear clean, standard laboratory gloves or use forceps when you handle them. Handling without gloves could contaminate the stones with oil, bacteria etc from skin.

Also, do not use powdered gloves, as these can introduce contaminants to the mobile phase.

Click [HERE](#) for ordering information and pictures of Sparging Stones.

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com