



Can amber or clear polypropylene vials be sterilized with autoclaving or ethylene oxide – FAQ

Polypropylene autosampler vials [9502S-PP-Clear](#) should not be autoclaved because they may warp during heating. However, you can use ethylene oxide gas to sterilize the vials if you do not exceed 120°C.

Click [HERE](#) for more information about polypropylene vials and pictures

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com