

Our filter / sparging stones have an internal thread of 5-40. These should be used with low pressure 1/8" or 1/16" ID polymeric tubing. The threads are actually small barbs that grab onto the tubing, ensuring a secure fit.

*NOTE: It may be important to use the internal threaded sparging stones instead of those with fittings to avoid the mixture of many different components in the solvent. This type of connection of the tubing and the stone should not come off in situ as can happen with the tubing/stem configuration.*

Click [HERE](#) for filter sparging stone ordering information and pictures

Click [HERE](#) for low pressure tubing ordering information and pictures

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)