
What are the maximum temperature and pressure specifications for the PEEK column hardware - FAQ

The PEEK columns can be used for short-term packing at pressures up to 500 bar (see [packing protocol](#)) and 300-350 bar under normal long-term operating conditions.

You can use the columns at temperatures up to 150-200 °C, but the maximum recommended operating pressure will be lower at elevated temperatures. In this temperature range, we recommend not exceeding 50-75 **bar**.

Also please note to use only solvents which are PEEK-compatible (*i.e. will not extract or soften PEEK*).

[PEEK Column Hardware Ordering Information](#)