



How does the shorter column supplied in the HSQ kit affect the data compared to the longer column from the PQ Kit - FAQ

The HSQ Kit™ comes with a 4.6 x 75mm column that is used in the qualification procedure.

The older version of the the HSQ Kit™ (called the PQ Kit™) comes with a 4.6 x 150mm column instead. This difference will not affect the qualification data results and the columns can be used **interchangeably**. The only difference for the end-user is that the HPLC runs take much less time due to the shorter column length when using the HSQ Kit™.

Hence, this a great advantage over the older model.

Click [HERE](#) for HSQ & PQ kits ordering information and pictures.

Printed from the Chrom Resource Center

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 06-16-2024