



What type of fitting is best to use with the Cogent HPLC and LCMS columns - FAQ

We generally recommend the Direct Adaptive Di-Ad™ column connectors instead of a fitting that can be re-swaged each time one of our columns is removed (*both standard 4um and 2.0™ phases*) to enable a clean connection when changing HPLC columns. The Di-Ad™ allows proper and easy installation of Cogent columns as well as different columns without the hassles of swaged ferrules.

We do not recommend Di-Ad™ or PEEK fittings for UHPLC.

Click [HERE](#) for Di-Ad™ column connector information and pictures.