MICROS

How do I calculate plug length in Capillary Electrophoresis?

One possible way is to use the "Poiseuille equation" to calculate the plug length:

 $V = (\Delta P d^4 \pi t)/(128 \eta L)$

where V is volume of the injected plug, ΔP is pressure drop down the length of the capillary (pascals), d is the inside diameter (m), η is the fluid viscosity (Pascal-sec), *t* is time the pressure is applied (sec), and *L* is the length of the capillary (m). Here you solve for V to get the injected plug volume. From there, you simply use the cylinder volume equation to convert volume to length of the plug:

 $V = \pi r^2 L_{plug}$

 $L_{plug} = (V / (\pi r^2))$

 $L_{plug} = (V / (\pi (d/2)^2))0$



where r is the inside radius and L_{plug} is the length of the plug.

Printed from the Chrom Resource Center Copyright 2024, All Rights Apply **MicroSolv Technology Corporation** 9158 Industrial Blvd. NE, Leland, NC 28451 tel. (732) 380-8900, fax (910) 769-9435 Email: customers@mtc-usa.com Website: www.mtc-usa.com