

When you insert the guard column cartridge and close the holder, you may notice that the holder does not appear to close all the way.

This is normal and within the design specifications of the unit. With the cartridge placed in the holder, simply give the holder a quarter turn with a wrench and this should suffice to provide adequate sealing of the guard column.



Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com