

## Can the U-2D Micro-Sample management system base racks and inserts be centrifuged - FAQ

Yes, with the rack nested into the Base, our U-2D™ will function similarly to a typical 96-well micro titer plate system. Currently, we have confirmed normal operation up to 2,500 x g.

Click [HERE](#) for U-2D Micro-Sample Management System ordering information and pictures



**U-2D™ 96 Well Plates**  
MICROSOLV

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 07-22-2024