



Why is it not recommended to use FEP tubing above 80°C when the melting point is much higher - FAQ

FEP polymer has a melting point of 260°C and a general recommended maximum operating temperature during usage of around 200°C.

If pressure is applied in the tubing, however, the maximum temperature should be lowered accordingly. The maximum temperature during usage should be 80°C under these conditions. If no pressure is applied, the tubing can be used up to 200°C.

Click [HERE](#) for low pressure HPLC Tubing ordering information and pictures.

Printed from the Chrom Resource Center

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 07-22-2024