



How do I clean the PQ or HSQ kit HPLC column – FAQ

Since the column should only be used for test purposes, it should never become contaminated, and should require only a minimal 10-15 minute equilibration prior to starting the sequence. If it ever requires cleaning, flush for 15-30 minutes with 50:50 acetonitrile : water.”

Click [HERE](#) for PQ & HSQ Kit ordering information and pictures.

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com