


NMR structure confirmation data for Cogent UDC-cholesterol phases – Tech Information

The following solid state CP-MAS ^{13}C NMR spectrum was recorded for a Cogent UDC-Cholesterol™ stationary phase in order to elucidate the structure. The assignments for the different carbon groups are shown below. The spectrum highlights the unique bonding of these phases, which is not possible with conventional silica based stationary phases.

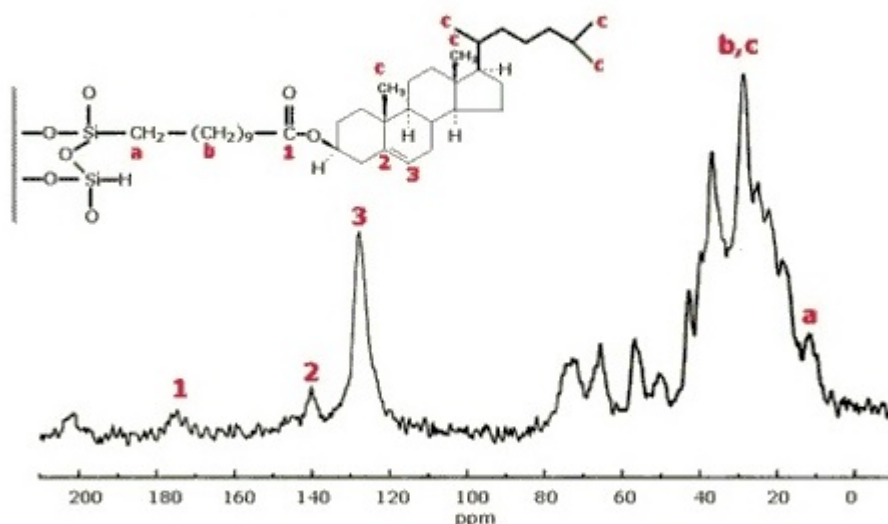

UDC-Cholesterol

Click [HERE](#) for Cogent UDC-cholesterol HPLC column ordering information



Printed from the Chrom Resource Center
MicroSolv Technology Corporation
9158 Industrial Blvd. NE, Leland, NC 28451
tel. (732) 380-8900, fax (910) 769-9435
Email: customers@mtc-usa.com
Website: www.mtc-usa.com
Date: 07-22-2024