

**You can distinguish U-2D™ inserts (used vs. unused) by the following ways:**

- A. Used samples can be indicated on the silicone mat/film using a marker.
- B. Alternatively, the film or mat can be cut with scissors to an appropriate size to protect unused inserts or to help keep track of used inserts.
- C. Users can selectively remove unused (or used) inserts as needed. Inserts should be carefully handled/stored to prevent contamination prior to use or transfer.

Click [HERE](#) for U-2D™ Micro-Sample Management System Ordering Information

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)