

Better methods than HILIC in HPLC and LCMS – Tips & Suggestions

Over the years, many chromatographers have been given the task to develop HPLC and LCMS methods for highly polar compounds. Those who have chosen HILIC columns have reported gradient methods can be problematic causing frustration, excessive method development time and additional costs. Mostly the issues are the “downtime” between runs that HILIC and all ordinary silica columns struggle with including equilibration of the on-column, silica water layer.

The permanent water layer or “hydration shell” causes difficulty maintaining precision and reproducibility especially after several runs or when laboratory conditions such as temperature changes. It’s also reported that the HILIC column lifetime can be short and the columns can fail without notice during over night runs.

High salt concentrations in the mobile phase (up to 100 mM) is often needed in HILIC to get proper retention times which can be a problem for LCMS instruments and preparative chromatography.

ANP or Aqueous Normal Phase HPLC is similar to HILIC in that polar compounds are retained and separated using high organic content mobile phases as a starting point and running an inverse gradient (high organic to high aqueous). A difference in that Cogent TYPE-C HPLC columns do not retain a water layer (see images below), does not have more than 5% residual silanols and utilizes direct silicon carbon bonds for ligands, the problems most often associated with HILIC are eliminated. Low salt concentrations are used in ANP (ie 15mM or less) and no buffering is needed thus avoiding all the problems of high salt concentrations in LCMS and preparative chromatography.

Some Benefits to Chromatographers using ANP v. HILIC:

- Faster Equilibration between runs (3-5 column volumes).
- Gain more Throughput per instrument, save on Cost of Solvents.
- Precision is Extraordinary run to run, day to day.
- More Confidence in your data, better for GMP and regulated laboratories.
- Column Lifetime increased over HILIC columns.
- Fewer failures during runs, lower cost, Budget Saver.
- Lower Salt needed for retention.
- Low Salt will not fowl the inlet of LCMS and offers faster recovery of Prep Samples.
- ANP can retain Polar & some non-Polar Compounds.
- Wider selectivity range, Fast Method Development.

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply


MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451


tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

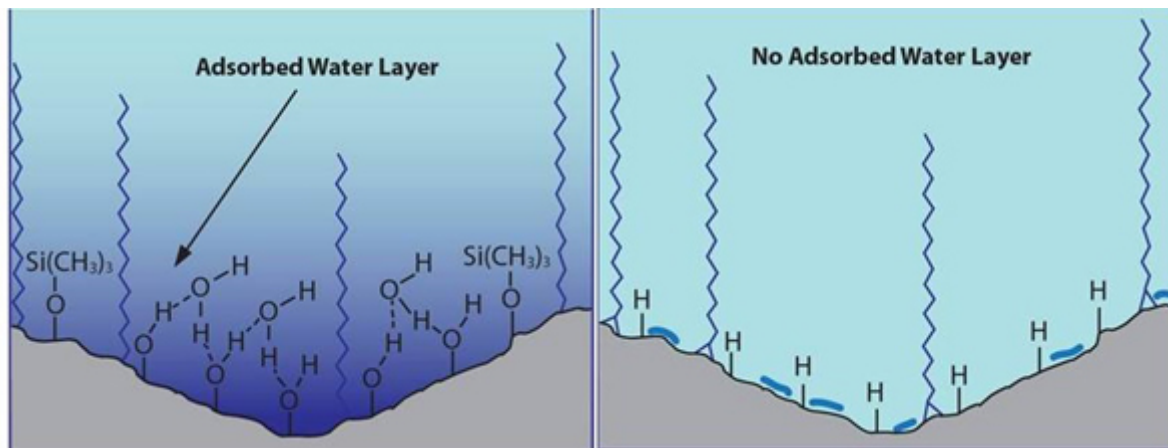

HILIC Phases have a Water Shell and Cogent TYPE-C Silica Does Not

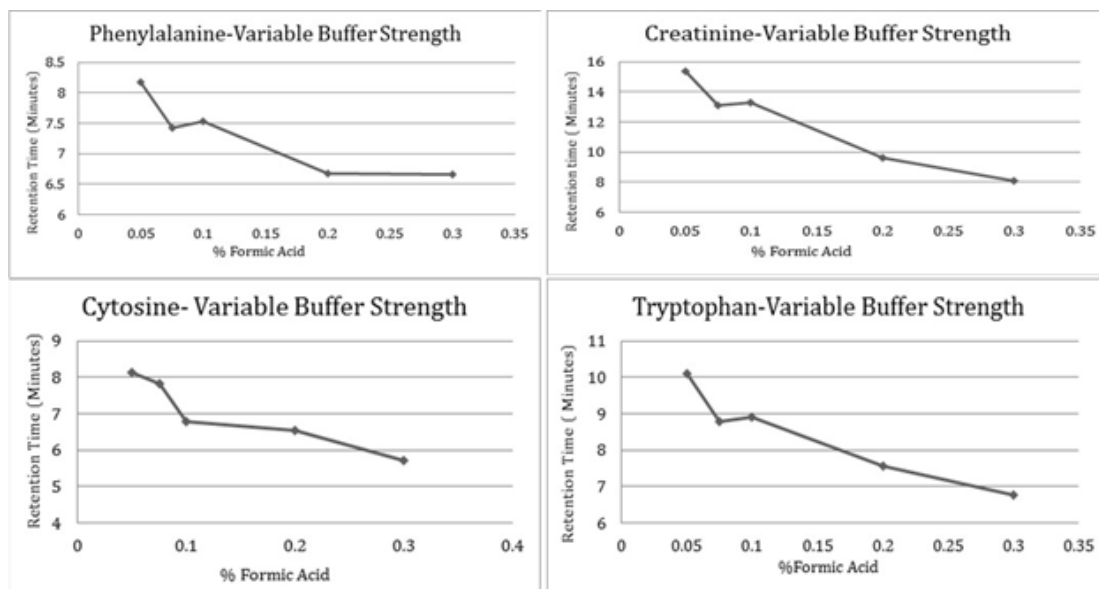

Effect of Salt Concentration on Retention in ANP Methods.

Notes:

1. Less than one Monolayer of Water on Cogent TYPE-C columns vs. 3-12 layers on HILIC materials leading to less variability in surface composition and hence greater Precision in analyte retention.
2. Rapid Equilibration of the stationary phase after gradients, typically in the range of 3-5 column volumes for Cogent TYPE-C Columns.
3. Negatively Charged Surface on Silica Hydride due to free, surface Hydroxide ions.
4. Difference in Selectivity and Enhanced Retention in ANP.
5. Ability of Cogent TYPE-C Columns to Function in Reversed-Phase HPLC Mode.
6. Excellent run to run and column to column Reproducibility with Cogent TYPE-C Columns.
7. No need for high Concentrations of Additives in the Mobile Phase; only 15mM or less required for most Hydrophilic Compound Analyses in ANP.

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451
tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

See also: [Main Differences between HILIC & Cogent TYPE-C Columns...](#)

See also: [Why ANP HPLC Methods are Better to use than HILIC.](#)

See also: [Efficiency of ANP v. HILIC Columns](#)

Reference below for details on this topic published in J. Phys. Chem. B.:

C. Kulsing, Y. Nolvachai, P.J. Marriott, R.I. Boysen, M.T. Matyska, J.J. Peseck, M.T.W. Hearn Insights into the Origin of the Separation Selectivity with Silica Hydride Adsorbents J. Phys. Chem. B. 2015 119 3063-3069



Attachment:

Very High Flow Rate, High Efficiency & Selectivity PDF [Download File](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com