



## Can autosampler vials be used in a centrifuge and what is the maximum centrifugal force that can be used – FAQ

**Autosampler vials are not designed to be centrifuged so we cannot support it with the exception of the MRQ™ & Max Recovery™ vials.**

Due to their thicker walls and “single piece” construction (AQ brand, RSA or RSA-Pro) they can withstand g forces up to 6,000 x g when used in a compatible centrifuge and adapters.

Click [HERE](#) for ordering information and pictures of the MRQ™ center draining vials.

Click [HERE](#) for ordering information and pictures of the Max Recovery™ vials.

Click [HERE](#) for ordering information and pictures of The Vial Centrifuge product page

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)