



Can a crimped cap be removed from Fused Insert vials without damaging the vial - FAQ

The answer depends on the type of de-capper that is used.

The type of de-capper with a rod that pushes downward onto the cap, as you squeeze the handles, into the vial will not be suitable, as this would damage the vial insert.

There is a type of de-dapper that grabs the cap from the sides, which resembles pliers and it quite dissimilar in appearance to a crimping tool, that can be used to carefully remove the cap. This holds true for vials with both fused and non-fused inserts.

MicroSolv does not sell this type of product.

Printed from the Chrom Resource Center

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 07-22-2024