



Can a crimped cap be removed from Fused Insert vials without damaging the vial - FAQ

The answer depends on the type of de-capper that is used.

The type of de-capper with a rod that pushes downward onto the cap, as you squeeze the handles, into the vial will not be suitable, as this would damage the vial insert.

There is a type of de-dapper that grabs the cap from the sides, which resembles pliers and is quite dissimilar in appearance to a crimping tool, that can be used to carefully remove the cap. This holds true for vials with both fused and non-fused inserts.

MicroSolv does not sell this type of product.

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com