



What is the main difference between ordinary and LCMS compatible polypropylene vials – FAQ

The differences between item **9532S-OPV** and **9532S-MS** is that item 9532S-MS is made from a special type of polypropylene and we do not use “mold release agents” in the molding process making it less likely to have any extractables or leachables under any circumstance.

Acetonitrile is likely to remove existing mold release agents which can show up in chromatograms as spurious peaks.

Printed from the Chrom Resource Center

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com

Date: 07-03-2024