





## Glycosidic Analyzed with LCMS – AppNote

### Cyanidine-3-Hexoside, Luteolin 7-O-Glucoside & Malvidin Dihexoside


Click [HERE](#) for Column Ordering Information.

Observed in Mesquite Flour extracts, these compounds are Glycosidic Flavones and Anthocyanins. The Glycoside Moieties impart polar characteristics to the analytes, which makes them amenable to Retention by the Aqueous Normal Phase (ANP).

In this Application Note, the Cogent Phenyl Hydride Column was used to obtain Chromatographic Separation. The same column can be used in the Reversed Phase mode as well, highlighting its versatility in the analysis of both hydrophilic and hydrophobic compounds.

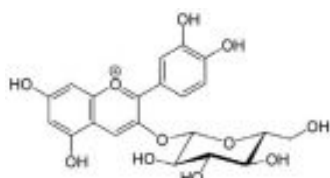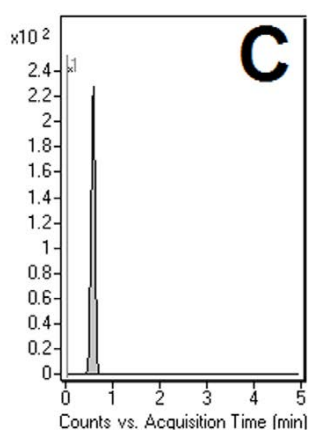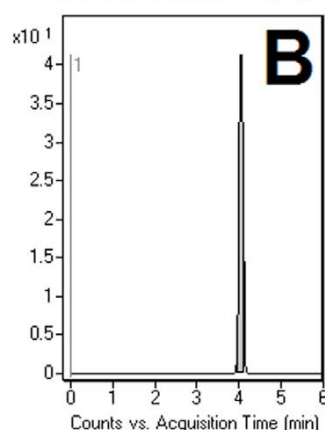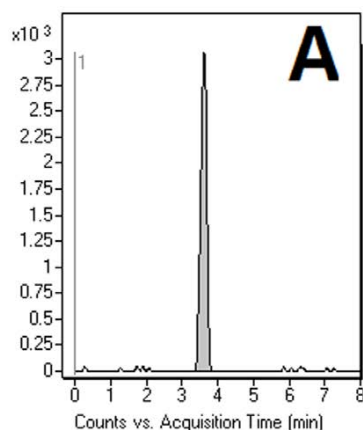

Cyanidine-3-hexoside

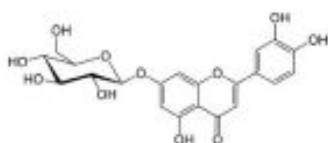

Luteolin 7-O-glucoside

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

**Peaks:**

A. Cyanidine-3-Hexoside ( $m/z = 450.116 [M + H]^+$ )

B. Luteolin 7-O-Glucoside ( $m/z = 331.082 [M + H]^+$ )

C. Malvidin Dihexoside ( $m/z = 449.108 [M + H]^+$ )

## Method Conditions:

**Column:** Cogent Phenyl Hydride™, 4 $\mu$ m, 100A

**Catalog No.:** 69020-05P-2

**Dimensions:** 2.1 x 50 mm

### Solvents:

A: DI Water / 0.1% Formic Acid (v/v)

B: Acetonitrile / 0.1% Formic Acid (v/v)

### Gradient:

Time (Minutes)	%B
0	90
10	50
12	50
18	20
20	20
21	90

**Injection Volume:** 1 $\mu$ L

**Flow Rate:** 0.4mL / minute

**Detection:** ESI - POS - Agilent 6210 MSD TOF Mass Spectrometer

**Sample Preparation:** Mesquite flour extracted in DI Water and 70:10 Methanol / DI Water. Extracts were centrifuged and the supernatant was collected.

**Note:** Mesquite Flour may have several advantages over wheat-based Flour for certain types of people. Firstly, it is gluten-free, which may be pertinent to those with celiac disease or gluten sensitivity. For diabetics, its low Glycemic index makes it a good source of Carbohydrates that that can help keep blood sugar levels stable. It is also high in protein and dietary fiber.



### Attachment

**No 370 Glycosidic Constituents of Mesquite Flour.pdf** 0.2 Mb [Download File](#)

Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)