

**Methanol is generally not recommended for most ANP applications** however a few scientists have seen success with it.

Instead, it is best to use polar aprotic solvents such as acetonitrile. If using ELSD or LCMS, acetone can also be a good alternative to acetonitrile as well. However, because of its low **UV cutoff**, it is generally not appropriate for analyses performed with UV detection.



Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)