


## Suggestions on how to separate capsaicinoids in hot sauce by HPLC – Tips & Suggestions

If you have a hot sauce extract containing nordihydrocapsaicin, capsaicin, dihydrocapsaicin, and homodihydrocapsaicin, the Cogent Bidentate C18 2.0™ HPLC column is a good choice for sufficient separation when the compounds are analyzed by LC-MS; look for the  $[M+H]^+$  ions.

The following gradient should be a good starting point:



You may get EICs which look like this:

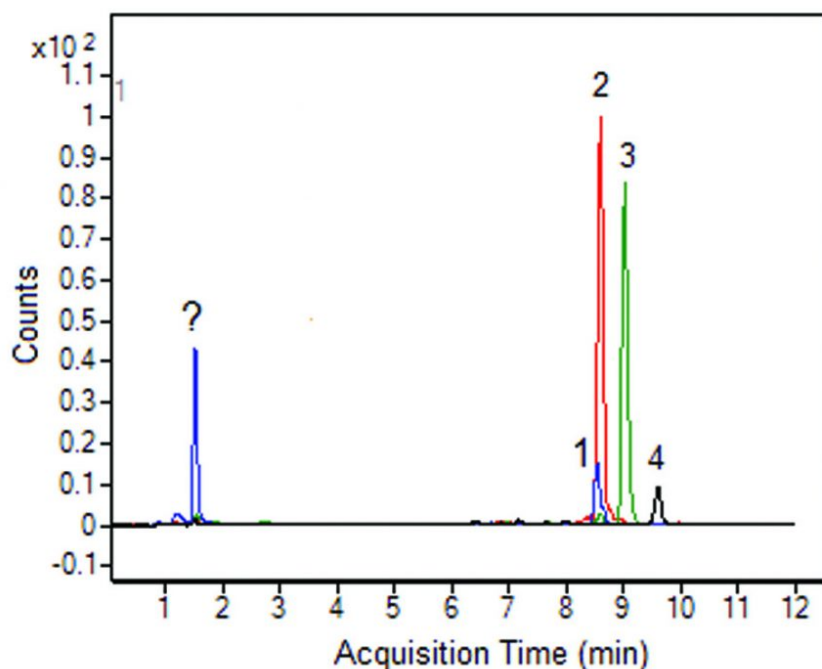

1. nordihydrocapsaicin, 294.2064 m/z  $[M+H]^+$

2. capsaicin, 306.2064 m/z  $[M+H]^+$

3. dihydrocapsaicin, 308.2220 m/z  $[M+H]^+$

4. homodihydrocapsaicin, 322.2377 m/z  $[M+H]^+$

For peaks 1 and 2, you will be able to obtain separate peaks for each compound in the EICs so there will not be an issue of interference from co-elution.

***In terms of sample preparation, try using a Soxhlet extraction procedure.***

[Click HERE for Cogent Bidentate C18 HPLC Column Ordering Information.](#)

# MICROSOLV

COGENT  
HPLC Columns™

Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 05-17-2024