

Since Cogent TYPE-C columns can be used in Reversed Phase or Normal Phase what do you suggest to switch these solvent systems which are not miscible - FAQ

Suggested Procedure:

A - Changing from Reversed Phase to Normal Phase HPLC; pump 100% methanol for 15 minutes at 1 mL / minute flow rate, followed by 15 minutes 100% methylene chloride. The column is ready to be equilibrated with mobile phase for NP HPLC.

B - Changing from Normal Phase to Reversed Phase HPLC; pump 100% methylene chloride for 15 minutes at 1 mL / minute flow rate, followed by 15 minutes 100% methanol. The column is ready to be equilibrated with mobile phase for RP-HPLC.

A simple 30 minute procedure allows switching from one mode to another.



Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

MicroSolv Technology Corporation

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: customers@mtc-usa.com

Website: www.mtc-usa.com