



## Change the type of methods used from DAD to VWD in the HSQ software – How To

1. To change the detector methods used from **DAD** to **VWD**, please remove data in **“Found Wavelengths from Scanning Detectors”** (~column 181) and reenter the values to the columns directly below, named: **“Found Wavelengths for Manual Determination”** (~column 185)
2. Please note, the values in **“Found Wavelengths from Scanning Detectors”** must be left blank with no data entered, otherwise it will print the final certificate with **DAD methods used**.



For more information, please read more [here](#).



Printed from the Chrom Resource Center

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 07-03-2024