

## Change the type of methods used from DAD to VWD in the HSQ software - How To

1. To change the detector methods used from **DAD** to **VWD**, please remove data in **“Found Wavelengths from Scanning Detectors”** (~column 181) and reenter the values to the columns directly below, named: **“Found Wavelengths for Manual Determination”** (~column 185)
2. Please note, the values in **“Found Wavelengths from Scanning Detectors”** must be left blank with no data entered, otherwise it will print the final certificate with **DAD methods used**.

**Data Entry for Detector Wavelength Assembly**

**Primary Standard (Minimum 0.05%)** Enter Retention/ Wavelength Data

Select up to 8 Retention/Wavelengths for Calibration from the following drug class lists, or specify the 8 official ICH bands

1.1	200 nm	1.2	210 nm	1.3	220 nm	1.4	230 nm	1.5	240 nm
-----	--------	-----	--------	-----	--------	-----	--------	-----	--------

Found Wavelengths for Scanning Detectors (see blank if Manual method is used)

1.1	200 nm	1.2	210 nm	1.3	220 nm	1.4	230 nm	1.5	240 nm
-----	--------	-----	--------	-----	--------	-----	--------	-----	--------

Found Wavelengths for Manual Determination

1.1	200 nm	1.2	210 nm	1.3	220 nm	1.4	230 nm	1.5	240 nm
-----	--------	-----	--------	-----	--------	-----	--------	-----	--------

**Secondary Standard (Optional)** Enter Calibration Data

Standard Calibration Wavelengths

1.1	200 nm	1.2	210 nm
-----	--------	-----	--------

Found Wavelengths for Scanning Detectors (see blank if Manual method is used)

1.1	200 nm	1.2	210 nm
-----	--------	-----	--------

Found Wavelengths for Manual Determination

1.1	200 nm	1.2	210 nm
-----	--------	-----	--------

Plumbridge

For more information, please read more [here](#).



Printed from the Chrom Resource Center

Copyright 2024, All Rights Apply

**MicroSolv Technology Corporation**

9158 Industrial Blvd. NE, Leland, NC 28451

tel. (732) 380-8900, fax (910) 769-9435

Email: [customers@mtc-usa.com](mailto:customers@mtc-usa.com)

Website: [www.mtc-usa.com](http://www.mtc-usa.com)

Date: 09-18-2024